

Polyclonal Production Protocol

Rabbit Protein Schedule of Events, 118 Day

Standard Schedule of Events

Day:	Procedure:
0	SPF Rabbit Pre-bleed (Avg. 5 ml serum) 1° SC: 250 micrograms with FCA
21	Boost SC: 125 micrograms with FIA
31	Test Bleed (Avg. 5 ml serum)
32-38	ELISA Titer Assay of Bleed *
42	Boost SC: 125 micrograms with FIA
52	Test Bleed (Avg. 5 ml serum)
63	Boost SC: 125 micrograms with FIA
73	Production Bleed (Avg. 20 ml serum)
84	Boost SC: 125 micrograms with FIA
94	Production Bleed (Avg. 20 ml serum)
95-101	ELISA Titer Assay of Bleed *
105	Boost SC: 125 micrograms with FIA
115	Production Bleed (Avg. 20 ml serum)
118	Terminal bleed (Avg. 50 ml serum)

Total material required for injection for the above program is 875 micrograms/animal.

If we are conjugating protein for you we assume a 20% loss of protein during conjugation.

Total amount of serum expected from a rabbit on the above schedule of events is approximately 120 ml. If sending gel format, please indicate amount of micrograms per vial or if indicating amount per gel slice, please indicate the number of gel slices.

Please Note:

- It is recommended that a minimum of two rabbits be used due to biological variations.
- Immunogen is emulsified in Freund's Complete Adjuvant (FCA) for initial injections. Freund's Incomplete Adjuvant (FIA) is used for all subsequent injections (boosts). Projects follow a three-week cycle of boosts. Test bleeds are taken approximately 10 days after the boosts.

***Optional ELISA:** We suggest testing the first bleed of the project to evaluate the effect of antigen on the animal. Information gained from an ELISA at this point will provide an opportunity for you to make adjustments to your immunization schedule. Additional ELISA's performed later in the project will give you valuable information that can be used to determine which sera samples to include in your research or which bleeds you would like to have purified.